



HandyStep® S

The uncomplicated repetitive pipette for complicated liquids

BRAND. For lab. For

- + High viscosity, high density, high vapor pressure – no problem, thanks to the positive displacement principle
- + Compatible with BRAND PD-Tips and other positive displacement tips
- + Robust, stable and functional

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Россия +7(495)268-04-70

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Казахстан +7(7172)727-132

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Киргизия +996(312)96-26-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93



Get to know the HandyStep® S

The HandyStep® S repetitive pipette combines the application advantages of the positive displacement principle with easy operation, robust design and an attractive price-to-performance ratio.

Thanks to the positive displacement principle, the HandyStep® S is often the easy solution for liquids that are difficult to pipette. This includes highly viscous or wetting media, high density media, foaming media such as detergents or highly volatile liquids.

Its balanced weight distribution and the practical ergonomic positioning of operating elements makes work more efficient, even when pipetting long series. The ergonomic finger rest

relieves strain on the hand, the positioning of the dispensing lever is easy on the thumb, and the stroke can be set with one hand, even while wearing gloves. Tips can be ejected with the press of a button, without having to touch the tips, making tip ejection contamination-free.

The HandyStep® S repetitive pipette works with the original BRAND PD-Tips II. Tips from many other manufacturers can also be used, providing maximum flexibility.

The simple, robust design and purely mechanical, battery-free operation make the HandyStep® S a reliable lab partner. Discover the application possibilities of the HandyStep® S.

HandyStep® S




- + Volume range from 2 µl to 5 ml
- + Adjustable to 59 volumes and up to 49 dispensing steps
- + Weighs just 108 g
- + Easy tip connection – simply insert the PD-Tip

At a glance: Advantages of the HandyStep® S

The HandyStep® S is a robust repetitive pipette that uses the positive displacement principle. With the PD-Tips, the HandyStep® S covers a wide volume range from 2 µl to 5 ml

and can be used in many different applications, including diagnostics, molecular biology and environmental analysis.

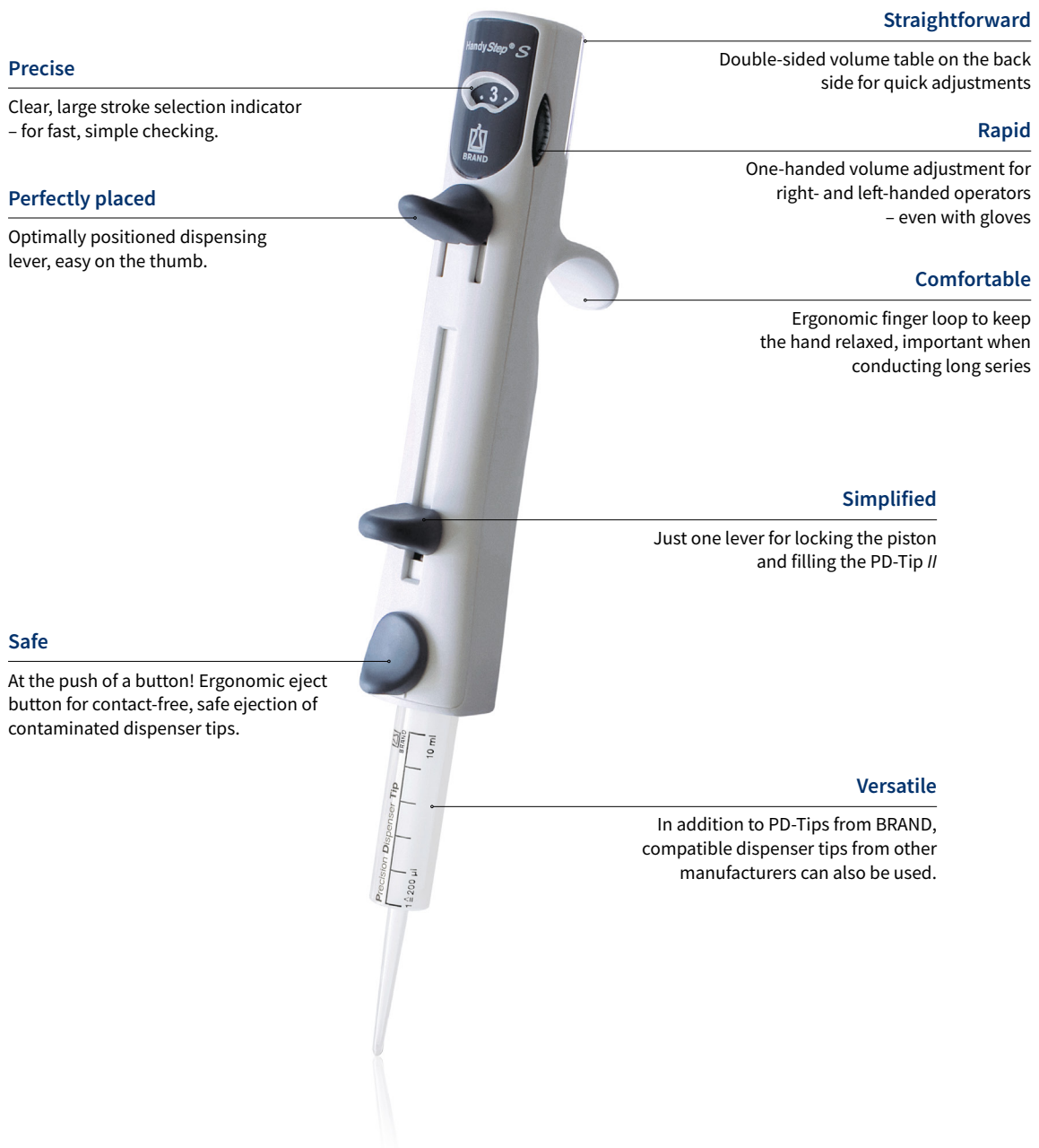

Volume table HandyStep® S with PD-Tips //

Offers 59 different partial volumes with different numbers of dispensing steps, depending on the PD-Tip size and stroke setting number used.

Setting	Tip size (ml)										Steps
	0.1	0.5	1	1.25	2.5	5	10	12.5	25	50	
1	2	10	20	25	50	100	200	250	500	1000	49
1.5	3	15	30	37.5	75	150	300	375	750	1500	32
2	4	20	40	50	100	200	400	500	1000	2000	24
2.5	5	25	50	62.5	125	250	500	625	1250	2500	19
3	6	30	60	75	150	300	600	750	1500	3000	15
3.5	7	35	70	87.5	175	350	700	875	1750	3500	13
4	8	40	80	100	200	400	800	1000	2000	4000	11
4.5	9	45	90	112.5	225	450	900	1125	2250	4500	10
5	10	50	100	125	250	500	1000	1250	2500	5000	9
Volume (µl)											



Repetitive dispensing



Dispensing small volumes



The rack mount also fits the Transferpette® S bench-top rack.



Fast volume selection with the volume table on the back side

Ordering information

HandyStep® S

repetitive pipette, DE-M marked


Items supplied:

performance certificate, operating manual, shelf/rack mount.



	Pack of	Cat. No.
	1	705110

Accessories for simple and efficient work

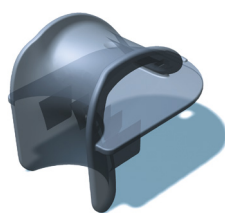

Shelf/rack mount

for 1 HandyStep® S repetitive pipette

	Pack of	Cat. No.
	1	705130



Wall mount

for 1 HandyStep® S repetitive pipette

	Pack of	Cat. No.
	1	705129

PD-Tips //

PD-Tips //, Precision dispenser tips with type code

Capacity ml	Pack of	non-sterile Cat. No.	BIO-CERT® LIQUID HANDLING STERILE Cat. No.
0.1	100	705700	705730
0.5	100	705702	705732
1	100	705704	705734
1.25	100	705706	705736
2.5	100	705708	705738
5	100	705710	705740
10	100	705712	705742
12.5	100	705714	705744
25	50 + 1 adapter	705716	705746*
50	25 + 1 adapter	705718	705748
Set	20 pieces per size: 0.5 1 1.25 2.5 5 10 12.5 ml	705720	



Adapter

for 25 and 50 ml PD-Tips //, PP,
autoclavable

Description	Pack of	Cat. No.
non-sterile	10	702398
BIO-CERT® LIQUID HANDLING STERILE	5	702399

* BIO-CERT®: pack of 25 + 1 adapter



Accuracy Table

The measurements were carried out with BRAND PD-Tips //. Distilled water was used as the test liquid. The nominal volume is the maximum volume printed on the PD-Tip. The tolerances specified in ISO 8655 are not exceeded.

PD-Tip //	Volume range	Accuracy A $\leq \pm \%$			Coefficient of variation CV $\leq \%$		
		Stroke setting $\triangleq$ % of nominal volume1 $\triangleq$			Stroke setting $\triangleq$ % of nominal volume1 $\triangleq$		
		1 $\triangleq$ 2%	3 $\triangleq$ 6%	5 $\triangleq$ 10%	1 $\triangleq$ 2%	3 $\triangleq$ 6%	5 $\triangleq$ 10%
0.1 ml	2 μ l - 10 μ l	4.0	2.4	1.6	6.0	3.0	2.0
0.5 ml	10 μ l - 50 μ l	2.5	1.5	1.0	2.5	1.5	1.0
1 ml	20 μ l - 100 μ l	2.5	1.5	1.0	2.0	1.2	0.8
1.25 ml	25 μ l - 125 μ l	2.5	1.4	0.9	2.0	1.1	0.7
2.5 ml	50 μ l - 250 μ l	1.8	1.1	0.7	1.5	0.9	0.6
5 ml	100 μ l - 500 μ l	1.8	1.1	0.7	1.5	0.9	0.7
10 ml	200 μ l - 1000 μ l	1.8	1.1	0.7	2.0	1.2	0.8
12.5 ml	250 μ l - 1250 μ l	1.8	1.1	0.8	3.2	2.0	1.4
25 ml	500 μ l - 2500 μ l	1.5	0.9	0.6	3.0	1.5	1.0
50 ml	1000 μ l - 5000 μ l	1.5	0.8	0.5	5.0	1.8	1.2

Алматы (7273)495-231	Иваново (4932)77-34-06	Магнитогорск (3519)55-03-13	Ростов-на-Дону (863)308-18-15	Тольятти (8482)63-91-07
Ангарск (3955)60-70-56	Ижевск (3412)26-03-58	Москва (495)268-04-70	Рязань (4912)46-61-64	Томск (3822)98-41-53
Архангельск (8182)63-90-72	Иркутск (395)279-98-46	Мурманск (8152)59-64-93	Самара (846)206-03-16	Тула (4872)33-79-87
Астрахань (8512)99-46-04	Казань (843)206-01-48	Набережные Челны (8552)20-53-41	Саранск (8342)22-96-24	Тюмень (3452)66-21-18
Барнаул (3852)73-04-60	Калининград (4012)72-03-81	Нижний Новгород (831)429-08-12	Санкт-Петербург (812)309-46-40	Ульяновск (8422)24-23-59
Белгород (4722)40-23-64	Калуга (4842)92-23-67	Новокузнецк (3843)20-46-81	Саратов (845)249-38-78	Улан-Удэ (3012)59-97-51
Благовещенск (4162)22-76-07	Кемерово (3842)65-04-62	Ноябрьск (3496)41-32-12	Севастополь (8692)22-31-93	Уфа (347)229-48-12
Брянск (4832)59-03-52	Киров (8332)68-02-04	Новосибирск (383)227-86-73	Симферополь (3652)67-13-56	Хабаровск (4212)92-98-04
Владивосток (423)249-28-31	Коломна (4966)23-41-49	Омск (3812)21-46-40	Смоленск (4812)29-41-54	Чебоксары (8352)28-53-07
Владикавказ (8672)28-90-48	Кострома (4942)77-07-48	Орел (4862)44-53-42	Сочи (862)225-72-31	Челябинск (351)202-03-61
Владимир (4922)49-43-18	Краснодар (861)203-40-90	Оренбург (3532)37-68-04	Ставрополь (8652)20-65-13	Череповец (8202)49-02-64
Волгоград (844)278-03-48	Красноярск (391)204-63-61	Пенза (8412)22-31-16	Сургут (3462)77-98-35	Чита (3022)38-34-83
Вологда (8172)26-41-59	Курск (4712)77-13-04	Петрозаводск (8142)55-98-37	Сыктывкар (8212)25-95-17	Якутск (4112)23-90-97
Воронеж (473)204-51-73	Курган (3522)50-90-47	Псков (8112)59-10-37	Тамбов (4752)50-40-97	Ярославль (4852)69-52-93
Екатеринбург (343)384-55-89	Липецк (4742)52-20-81	Пермь (342)205-81-47	Тверь (4822)63-31-35	
Россия +7(495)268-04-70	Казахстан +7(7172)727-132	Киргизия +996(312)96-26-47		

bdf@nt-rt.ru || <https://brand.nt-rt.ru/>